

INFORMATION ON

SETTING UP

THE SMARTRACK®



- Whenever using the SmartRack®, make sure that the SmartRack® or SmartRack® mini is placed on a stable and flat surface.
- Check the functionality of the SmartRack® and, if necessary, adjust the slot rails and module holders to an optimal position for your needs.

Assembly information for SmartRack® and SmartRack® mini by Better Basics Laborbedarf:


1. Begin by removing the SmartRack® or SmartRack® mini from its packaging and placing it on its feet. Make sure to place it on a flat surface so that it is stable and straight.






2. Note that each of the feet of the SmartRack® consists of two parts. The rear part of the feet can be removed by pulling it backwards, which allows the SmartRack® to be placed closer to a wall.



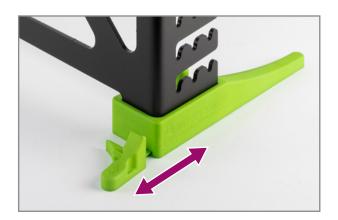

3. Place the drip tray on the two round positioning points on the feet.

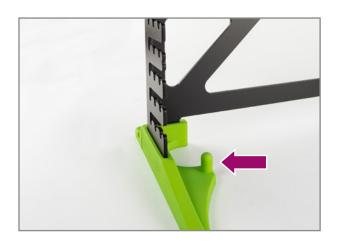




4. Hang the required number of slot rails (usually three to four slot rails) in the SmartRack® by holding them by the handles on the left and right side of the slot rail and move them to the desired position. There are three possible positions at each height. Please always change the position of the slot rails without the holder modules hooked in.





5. Insert the holder modules into the slot rails and then turn them from top to bottom until you feel a click. The holder modules are labelled with their respective names on the top. Make sure that each holder module is correctly positioned (clicked into place). A picture of the click mechanism is shown schematically here.



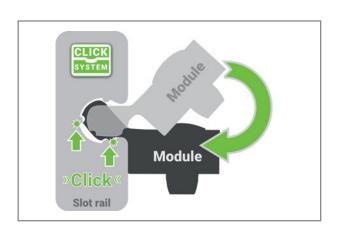



Once all the holder modules are positioned in the SmartRack®, you can hang the laboratory equipment. Hang the individual parts of your work equipment in the holder module provided.





7. Some holder modules have an additional, rotatable safety mechanism that protects the mounted laboratory glassware from falling out. To secure the equipment, turn the black locking ring until the ring closes the open front part of the holder module fully. To open the safety mechanism, proceed in reverse order.





8. With these steps, you have now successfully assembled your SmartRack® or SmartRack® mini and you are ready to store your equipment in a safe and orderly manner.

YOU ARE INTERESTED IN OUR PRODUCTS **OR YOU HAVE FURTHER QUESTIONS?**

Get in touch with us: contact@better-basics.de

You can also find more information and news about our product innovations at www.Better-Basics-Laborbedarf.de

Better Basics Laborbedarf GmbH

Loebtauer Str. 69, 01159 Dresden, Germany

- T: +49 (0) 178 9617 577 (Direct customers)
- +49 (0) 176 6233 8026 (Sales partners) T:
- E: contact@better-basics.de
- w: www.Better-Basics-Laborbedarf.de





