



# Product catalogue

## SmartRack®

Mobile Organisational Solutions for Laboratory Workstations



# Table of Contents

Your Advantage in Everyday Laboratory Work.....	3
SmartRack®   Mobile Organisational System.....	6
The Configuration.....	6
Frame size and type.....	7
Pre-configurations.....	9
Accessories.....	19
SmartCase.....	22
Customising.....	23
Item Overview.....	25
Technical Information.....	26

# Your Advantage in Everyday Laboratory Work

## Better Basics Laborbedarf – Your Advantage in Everyday Laboratory Work

Better Basics Laborbedarf optimises laboratories and their workspaces with innovative, award-winning organisational systems. We support our customers in creating the best possible working environment for their laboratory staff, helping to unlock the full potential of their laboratories. This approach ensures consistently high quality, efficiency, and process stability in the lab while providing a safe, ergonomic workplace with satisfied employees.

### Enhancing Employee Satisfaction

Our SmartRack® and SmartIntegrate product families offer a comprehensive system to optimise laboratory workflows. With over 100 modular solutions, they not only ensure order but also noticeably improve working conditions. Enhanced ergonomics, reduced movement paths, and clearly structured workstations increase both efficiency and employee satisfaction. Safe and organised environments reduce stress and promote job satisfaction. Furthermore, they demonstrably contribute to improved workplace safety and occupational health through thoughtful placement and ease of use.


### A System for Sustainable Process Improvement

Our product families are designed to standardise workflows and save time. The modular concept of SmartRack® and SmartIntegrate allows for flexible adaptations to any laboratory environment – from safety workbenches to large open-plan labs. These systems not only optimise space utilisation but also increase efficiency by an average of 30% and save up to 80% space. With a wide range of modules and accessories, they offer customised solutions for every need.

### Efficient, Safe, and Satisfying

For decision-makers and managers, SmartRack® and SmartIntegrate ensure process stability and economical resource use. At the same time, these systems provide innovation managers and organisational developers with the ideal basis for forward-looking workplace designs. For laboratory staff, this means fewer errors thanks to well-structured work areas, a more relaxed working environment, and often the ability to finish earlier due to significantly improved workflows.

**Discover how Better Basics Laborbedarf can transform your everyday laboratory work – for more efficient processes, satisfied employees, and excellent results.**

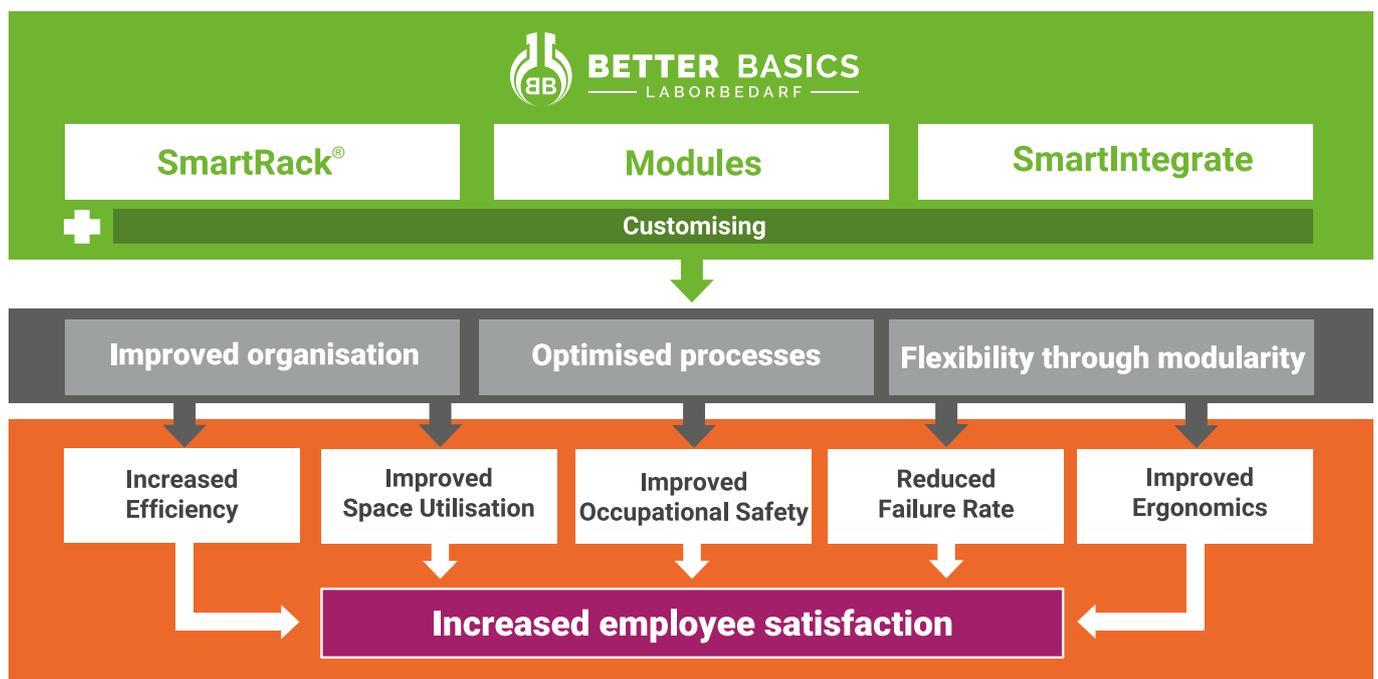

## The product overview of Better Basics

### ▶ SmartRack®, the mobile solution for: laboratory workbench, fume hood and safety cabinet



#### + Customised solutions

### ▶ SmartIntegrate, the integrative solutions for

#### I. Laboratory workbench



#### II. Fume hood



#### III. Safety cabinet



#### IV. Integration into laboratory equipment



#### + Customised solutions

### ▶ Modules

- Currently, more than 100 modules
- All modules in the Better Basics catalogue can be used universally in all SmartRack® and SmartIntegrate solutions
- Customised modules can be developed based on the specific customer needs, all customised modules can be used in all SmartRack® and SmartIntegrate solutions



#### + Customised solutions

▶ **SmartRack®**, the mobile solution for: laboratory workbench, fume hood and safety cabinet



✚ Customised solutions

## The SmartRack® – Efficiency and Safety Conveniently Combined.

The SmartRack® is the universal solution for laboratory organisation. With its modular and flexible structure, it optimises workspace utilisation and achieves up to 80% space savings. Whether in a laboratory fume hood, on the workbench, or in a biosafety cabinet – the SmartRack® adapts to your needs.

It supports efficient workflows, minimises sources of error, and enhances safety. Its pre-configured variants and customisable elements make the SmartRack® the ideal foundation for any laboratory environment.



Photo: A SmartRack® classic in a fume hood with tablet and laboratory equipment

# SmartRack® | The Configuration

<p><b>1. Selection of pre-configurations:</b> Basic Equipment, Biology, Chemistry, Pipetting Set, Consumables, etc.</p>			
<b>OR</b>			
<p><b>2. Customised configuration:</b> 2.1 Selection of frame size and frame type: SmartRack® primus (mini), SmartRack® classic (mini)</p>			
<p>2.2 Selection of individual modules: Holders and storage boxes as well as storage box inserts</p>			
<p>2.3 Selection of additional SmartRack® accessories (optional): If required further slot rail, drip tray or protective cover</p>			

## Example configuration of a personalised SmartRack® for Item no. 01704 and 01707, SmartRack® (left) plus modules (right)



Photo above: SmartRack® classic

Photos above: Selection of modules

**Each SmartRack® classic (Item number 01704) or SmartRack® classic mini (Item number 01707) is already delivered ex works with the following accessories**

- 1 x Pair of feet
- 1 x Drip tray
- 3 x Slot rails with handles
- 1 x Frame

# SmartRack® | Frame size and type

## SmartRack® classic



The SmartRack® classic was developed for use at many different laboratory workstations and especially for work in fume hoods. The SmartRack® classic mini is specially designed for work in safety cabinets. It consists of an aluminium frame with a black epoxy-polyester-based powder coating. Every third slot rail holder on both sides of the SmartRack® classic frame has a recess at the front to make it easier to align the slotted rails horizontally. The SmartRack® classic and the SmartRack® classic mini include the following components:

- 1 x frame incl. one pair of rear shortenable feet with tray holder
- 3 x slot rails with handles
- 1 x drip tray

Size comparison SmartRack® classic (left) and SmartRack® classic mini (right):



Item	Type	PU	Item no.	Price in EUR
SmartRack® classic	635 x 615 x 265 mm (WxHxD)	1	01704	699,00
SmartRack® classic mini	450 x 380 x 265 mm (WxHxD)	1	01707	499,00

The specified width of the SmartRack® and SmartRack® mini includes the frame, the slot rails with handles, as well as the feet.

If you need a SmartRack® with special size requirements or custom brackets, please contact us at [contact@better-basics.de](mailto:contact@better-basics.de).

## SmartRack® primus



The SmartRack® primus is the affordable entry-level version into the smart world of laboratory equipment from Better Basics. It was developed for use at many different laboratory workstations and in particular for work at the fume hood. The SmartRack® primus mini is specially designed for working on the safety cabinet. It consists of an uncoated, polished aluminium frame. Every third slot rail holder on both sides of the SmartRack® primus frame has a recess at the front to make it easier to align the slotted rails horizontally. The SmartRack® primus and the SmartRack® primus mini comprise the following components:

- 1 x frame incl. one pair of aluminium feet
- 2 x slot rails with handles

Size comparison SmartRack® primus (left) and SmartRack® primus mini (right):



Item	Type	PU	Item no.	Price in EUR
SmartRack® primus	635 x 615 x 265 mm (WxHxD)	1	01711	449,00
SmartRack® primus mini	450 x 380 x 265 mm (WxHxD)	1	01712	299,00

The specified width of the SmartRack® primus and SmartRack® primus mini includes the frame, the slot rails with handles, as well as the feet.

If you need a SmartRack® with special size requirements or custom brackets, please contact us at [contact@better-basics.de](mailto:contact@better-basics.de).

# SmartRack® | Pre-configurations

## SmartRack® mini pipetting set



The SmartRack® mini pipetting set has been specially designed for working with liquids in small quantities and includes the following components (item number of the individual components in brackets for better understanding):

9 components:

- 1 x SmartRack® classic mini (Item no. 01707)
- 1 x Holder for bin liner (Item no. 01202)
- 1 x Holder for pipettes - for 3 pipettes with hook (Item no. 01145)
- 2 x Holder for pipette tip boxes (Item no. 01147)
- 2 x Holder for laboratory screw top bottles (Item no. 01122)
- 2 x Holder for well plates (Item no. 01125)

Detailed views of the individual components included:



Item no. 01707



Item no. 01202



Item no. 01145



Item no. 01147



Item no. 01122



Item no. 01125

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de). There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Laboratory screw top bottles, pipettes and pipette tip boxes, waste bags and well plates are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		9	var.	968,92
Price advantage:				-69,92
SmartRack® mini pipetting set, bundle price:		1	01512	899,00

## SmartRack® basic equipment



The SmartRack® basic equipment is a basic set for organising basic accessories in everyday laboratory use and includes the following components (item number of the individual items in brackets for better understanding):

6 components:

- 1 x SmartRack® (Item no. 01704)
- 1 x Storage box (Item no. 01142)
- 1 x Holder for bin liner (Item no. 01202)
- 1 x Holder for glove / tissue boxes (Item no. 01143)
- 1 x Holder for wash bottles (Item no. 01146)
- 1 x Holder for kitchen roll (Item no. 01162)

Detailed views of the individual components included:



Item no. 01704



Item no. 01202



Item no. 01142



Item no. 01143



Item no. 01146



Item no. 01162

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de). There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Waste bags, tissue boxes, kitchen paper rolls and wash bottles are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		6	var.	1.078,95
Price advantage:				-79,95
SmartRack® basic equipment, bundle price:		1	01506	999,00

## SmartRack® starter set Biology



The SmartRack® starter set Biology has been specially designed for work in the field of biology and includes the following individual components (item number of the individual component in brackets for better understanding):

8 components:

- 1 x SmartRack® (Item no. 01704)
- 2 x Holder for laboratory screw top bottles (Item no. 01122)
- 2 x Storage box (Item no. 01142)
- 1 x Holder for pipettes - for 3 pipettes with hook (Item no. 01145)
- 1 x Holder for pipette tip boxes (Item no. 01147)
- 1 x Holder for microlitre syringes (Item no. 01155)

Detailed views of the individual components included:



Item no. 01704



Item no. 01122



Item no. 01142



Item no. 01145



Item no. 01147



Item no. 01155

### Note on the Item numbers provided:


The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de).

There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Laboratory screw top bottles, pipettes, pipette tip boxes and microlitre syringes are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		8	var.	1.203,93
Price advantage:				-104,93
SmartRack® starter set Biology, bundle price:		1	01507	1.099,00

## SmartRack® starter set Chemistry

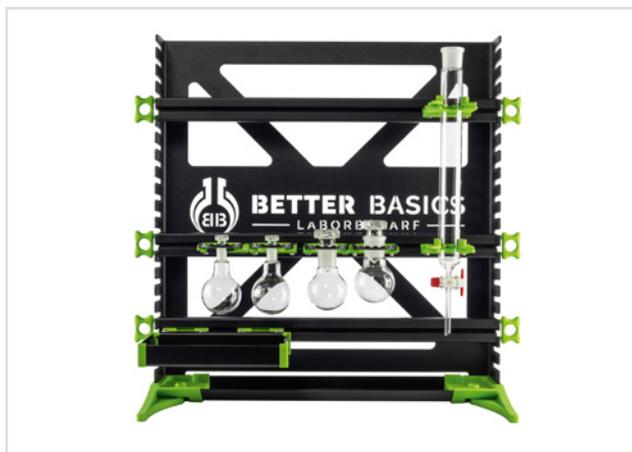

The SmartRack® starter set Chemistry has been specially designed for work in the field of chemistry and includes the following components (item number of the individual components in brackets for better understanding):

8 components:

- 1 x SmartRack® (Item no. 01704)
- 2 x Holder for glassware with ST14, plus closure (Item no. 01112)
- 2 x Holder for glassware with ST29, plus closure (Item no. 01115)
- 2 x Holder for chromatography column (Item no. 01107)
- 1 x Storage box (Item no. 01142)

Detailed views of the individual components included:



Item no. 01704



Item no. 01112



Item no. 01115



Item no. 01107



Item no. 01142

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de).

There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Chromatography columns and glassware are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		8	var.	1.178,93
Price advantage:				-79,93
SmartRack® starter set Chemistry, bundle price:		1	01508	1.099,00

## SmartRack® for consumables



The SmartRack® for consumables has been specially designed for the provision of consumables at the laboratory workplace and comprises the following components (item number of the individual components in brackets for better understanding):

12 components:

- 1 x SmartRack® (Item no. 01704)
- 1 x Dispenser for Parafilm® rolls (Item no. 01171)
- 1 x Dispenser for adhesive tapes (Item no. 01172)
- 3 x Holder for glove / tissue boxes (Item no. 01143)
- 1 x Holder for bin liner (Item no. 01202)
- 3 x Holder for wash bottles (Item no. 01146)
- 1 x Holder for kitchen roll (Item no. 01162)
- 1 x Slot rail with handles (Item no. 01601)



Detailed views of the individual components included:



Item no. 01704



Item no. 01171



Item no. 01172



Item no. 01143



Item no. 01202



Item no. 01146



Item no. 01162



Item no. 01601

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de).

There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Parafilm® rolls, adhesive tapes, glove/tissue boxes, waste bags, wash bottles and kitchen rolls are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		12	var.	1.673,89
Price advantage:				-174,89
SmartRack® for consumables, bundle price:		1	01509	1.499,00

## SmartRack® for Life Science



The SmartRack® for Life Science has been specially designed for work in the fields of biology, biomedicine, biotechnology as well as special fields of biology and for handling liquids in small quantities and preparing and storing samples for later analysis. It includes the following components (item number of each component in brackets for better understanding):

16 components:

- 1 x SmartRack® (Item no. 01704)
- 1 x Holder for bin liner (Item no. 01202)
- 1 x Storage box (Item no. 01142)
- 1 x Storage box insert - for small accessories (Item No. 01148)
- 1 x Storage box insert - for microreaction vessels (Item no. 01130)
- 1 x Storage box insert - for centrifuge tubes, (Item no. 01102)
- 2 x Holder for laboratory screw top bottles (Item no. 01122)
- 2 x Holder for well plates (Item no. 01125)
- 1 x Holder for wash bottles (Item no. 01146)
- 1 x Holder for pipettes - for 3 pipettes with hook (Item no. 01145)
- 2 x Holder for pipette tip boxes (Item no. 01147)
- 1 x Dispenser for Parafilm® rolls (Item no. 01171)
- 1 x Slot rail with handles (Item no. 01601)



Detailed views of the individual components included:



Item no. 01704



Item no. 01202



Item no. 01142



Item no. 01148



Item no. 01122



Item no. 01125



Item no. 01146



Item no. 01145



Item no. 01147



Item no. 01171



Item no. 01130



Item no. 01102



Item no. 01601

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de).

There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Laboratory screw top bottles, small accessories, pipette tip boxes, pipettes, microreaction vessels, sample vials, well plates, waste bags, Parafilm® rolls and wash bottles are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		16	var.	1.713,85
Price advantage:				-164,85
SmartRack® for Life Science, bundle price:		1	01510	1.549,00

## SmartRack® for Chromatography



The SmartRack® for Chromatography has been specially designed for substance separation tasks by means of column chromatography and comprises the following components (item number of the individual components in brackets for better understanding):

17 components:

- 1 x SmartRack® (Item no. 01704)
- 1 x Storage box (Item no. 01142)
- 1 x Storage box insert – for small accessories (Item no. 01148)
- 1 x Storage box insert - for sample containers and test tubes, 12x Ø 10 - 20 mm (Item no. 01160)
- 1 x Storage box insert - for sample containers and test tubes, 20x Ø 3 - 10 mm (Item no. 01161)
- 6 x Holder for chromatography column (Item no. 01107)
- 1 x Holder for wash bottles (Item no. 01146)
- 1 x Holder for microliter syringes (Item no. 01155)
- 2 x Holder for laboratory screw top bottles (Item no. 01122)
- 1 x Holder for glove / tissue boxes (Item no. 01143)
- 1 x Slot rail with handles (Item no. 01601)



Detailed views of the individual components included:



Item no. 01704



Item no. 01142



Item no. 01107



Item no. 01146



Item no. 01155



Item no. 01122



Item no. 01148



Item no. 01160



Item no. 01161



Item no. 01143



Item no. 01601

**Note on the Item numbers provided:**

The modules listed here can be found in the Better Basics partial catalog "Modules". You can download this catalog at [www.download.better-basics-laborbedarf.de](http://www.download.better-basics-laborbedarf.de).

There, you will also find the partial catalog SmartIntegrate as well as numerous additional details about our products and solutions.

Chromatography columns, wash bottles, microlitre syringes, laboratory screw top bottles, small accessories, tissue boxes, sample containers and microreaction vessels as well as glassware and glass devices are not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Price of the individual modules:		17	var.	1.823,84
Price advantage:				-124,84
SmartRack® for Chromatography, bundle price:		1	01511	1.699,00

## SmartRack® | Accessories

### Drip tray



The drip tray prevents the spread of liquids. In case of contamination, the drip tray can be easily removed from the SmartRack® and cleaned.

Item	Type	PU	Item no.	Price in EUR
Drip tray	for SmartRack®	1	01603	99,99
Drip tray mini	for SmartRack® mini	1	01604	89,99


### Slot rail with handles



The slot rail can be used in many different positions in the SmartRack®. A larger number of slot rails offers more options for placing modules through an optimal vertical arrangement, providing more scope for customizing the structure of the SmartRack® and adapting the module arrangement to the individual needs of each user.

Item	Type	PU	Item no.	Price in EUR
Slot rail with handles	for SmartRack®	1	01601	74,99
Slot rail mini with handles	for SmartRack® mini	1	01602	54,99

## Back-to-back connectors for SmartRack®

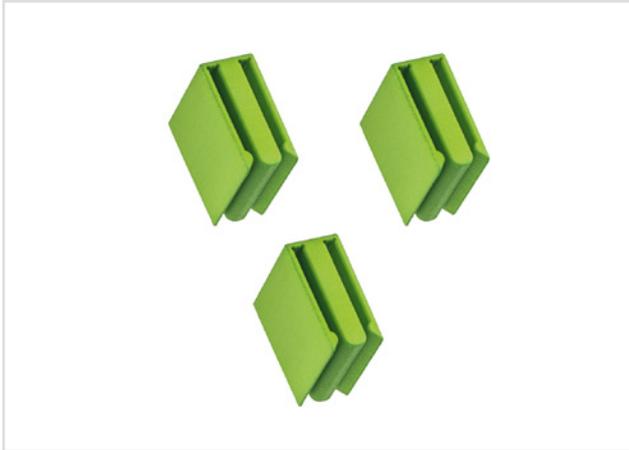

The back-to-back connectors for the SmartRack® enable two SmartRack® to be securely connected and thus used as a isolated solution. With the rear foot sections removed, two SmartRack® can be securely connected by simply attaching the connectors to the horizontal or diagonal braces of the frame rear panel. A set consists of three individual back-to-back connectors.

Item	Type	PU	Item no.	Price in EUR
Back-to-back connectors for SmartRack®		1	01614	79,99

## Protective cover for SmartRack®



Protective cover for SmartRack® with zipper\*

The protective hood made of fiber-reinforced polyethylene (mesh film), which was specially designed for the SmartRack®, reliably protects the SmartRack® against dirt and dust particles as well as providing splash protection against aggressive liquids. The dust and particle protection hood is available in two versions that do not differ in size. As a basic model and as a model with two zippers on the edges of the front. For regular removal of the dust and particle protection hood and for use in unfavorable height conditions, we recommend choosing the model with zippers.



Protective cover for SmartRack® without zipper\*



Protective cover for SmartRack® mini with zipper\*

\*) SmartRack® and equipment is not included in the scope of delivery.

Item	Type	PU	Item no.	Price in EUR
Protective cover for SmartRack®	with zipper	1	01609	139,99
Protective cover for SmartRack®	without zipper	1	01608	99,99
Protective cover for SmartRack® mini	with zipper	1	01613	89,99
Protective cover for SmartRack® mini	without zipper	1	01612	64,99

# SmartCase

## SmartCase



The SmartCase is designed to transport and store the SmartRack®, a modular organization system for laboratory workstations. Its aluminium case safely protects a SmartRack® inside and its contents from light and other environmental influences.

The SmartCase consists of a lower part with four castors into which a SmartRack® can be placed. The upper part of the SmartCase has a cushion on the inside, which serves to fix the SmartRack® in the SmartCase as soon as the SmartCase is closed by the tension locks. This prevents the SmartRack® from slipping or tipping over.

The SmartCase has a self-acting valve system which can be used to change the atmosphere inside the SmartCase by means of an external gas supply. This allows, for example, the temporary generation of a low-oxygen atmosphere in the SmartCase, which creates an additional safety level when transporting oxygen-sensitive, highly flammable substances or substances that spontaneously ignite in air.

For further information on the durability of the atmosphere exchange in the SmartCase, please contact us at [contact@better-basics.de](mailto:contact@better-basics.de).

The SmartCase includes the following components:

- A bottom section with four casters, four tension locks and an automatic gas outlet valve system
- A top part with a handle, four put-on aids and an automatic gas inlet valve system
- A connector for connecting gas hoses of various diameters to the SmartCase

*SmartRack® and other equipment are not included in the scope of delivery.*

Item	Type	PU	Item no.	Price in EUR
SmartCase		1	02101	949,00

# Customising

Better Basics Laborbedarf offers an unparalleled approach to laboratory workstation organisation with its customisable and modular solutions. Below is an overview of how our customisation services span both the SmartRack® and SmartIntegrate product lines:

## 1. Customisation of the SmartRack®

### For the SmartRack®

Our SmartRack® system is fully customisable to meet the specific requirements of your workspace. The dimensions of the SmartRack® can be individually adapted to the size of your equipment and workflows, ensuring seamless integration into your laboratory workstation. At the same time, the modular compatibility and reliability of your tailor-made SmartRack® are guaranteed through appropriately dimensioned slot rails. From compact configurations to larger frames, we enhance the organisation and efficiency of your laboratory.

## 2. Customisation of SmartIntegrate Products

### a) Bespoke SmartFrame for Fume Hoods

The SmartFrame optimises the organisation within fume hoods, offering a structured and safe arrangement for laboratory glassware, tools, and equipment. Custom-built SmartFrames provide a perfect fit for your specific workspace.

### b) Tailored SmartRail for Lab Benches

SmartRails bring adaptability and precision to your laboratory benches. Designed for both high-throughput workflows and compact setups, they ensure reduced clutter and enhanced efficiency. Each SmartRail is tailored to your preferred length and mounting specifications.

### c) Custom Solutions for Safety Cabinets

For safety cabinets, our custom solutions include integrated SmartRails, SmartFrames and SmartEdges designed to maximise space utilisation while supporting your workflows. Each component is built to ensure easy cleaning and adherence to safety regulations.

## 3. Custom Modules

### Tailored for any Laboratory Equipment

Using advanced 3D printing technology, we create modules and holders in any shape or size to fit your laboratory equipment. From pipette holders to mounts for specialised glassware, our bespoke designs cater to your unique requirements. Our rapid prototyping process ensures swift delivery of tailored solutions.

### Why Choose Better Basics for Customisation?

1. **Precision Craftsmanship:** Each component is tailored to your exact specifications, guaranteeing optimal performance and seamless integration.
2. **Fast Delivery:** Our in-house design and manufacturing teams ensure swift turnaround from concept to delivery.
3. **Transparent Pricing:** Costs for customisation are agreed upon in advance, providing clarity and control.
4. **Scalability:** Whether expanding an existing laboratory or building a new one, our products adapt to meet your growing needs.

### Transform Your Laboratory Workspace

With Better Basics Laborbedarf, your laboratory can benefit from customised SmartRack® or SmartIntegrate products.

Visit [customising.better-basics-laborbedarf.de](https://customising.better-basics-laborbedarf.de) or contact our team today to unlock the full potential of your laboratory space.

This concise yet comprehensive approach ensures your laboratory benefits from enhanced efficiency, organisation, and safety—all tailored to your specific needs.

## Examples for realised customisations

### Special frame sizes

#### SmartRack® frame - customisable in width as well as height



#### SmartRack® frame - extra wide

For a special requirement we manufactured an extra wide frame of the SmartRack®. Due to the width, this SmartRack® makes optimal use of the working space in our customer's laboratory.

### Special module sizes

#### Holder for multiple glassware



#### Double holder for glassware with ST10

For a project, a customer needed a double holder for glassware ST10. The module allows glassware with standard ground joint ST10 to be easily hung next to each other. The specially shaped recess prevents the glassware from falling out.

Of course, adaptations of this kind are also conceivable for other modules, e.g. storage boxes in individual sizes.

### Special laboratory equipment

#### Dispenser for cylindrical Petri dish multi-packs



For a globally operating biotechnology company, a dispenser for cylindrical Petri dish multi-packs was developed.

The dispenser not only enables easy removal of the packs but also provides a quick overview of the remaining stock.

# Item overview

## SmartRack®

Item	Type	Item no.	PU	Price in EUR	Page
SmartRack® classic	635 x 615 x 265 mm (WxHxD)	01704	1	699,00	7
SmartRack® classic mini	450 x 380 x 265 mm (WxHxD)	01707	1	499,00	7
SmartRack® for Chromatography		01511	1	1.699,00	17
SmartRack® for Life Science		01510	1	1.549,00	14
SmartRack® for consumables		01509	1	1.499,00	13
SmartRack® basic equipment		01506	1	999,00	10
SmartRack® mini pipetting set		01512	1	899,00	9
SmartRack® primus	635 x 615 x 265 mm (WxHxD)	01711	1	449,00	8
SmartRack® primus mini	450 x 380 x 265 mm (WxHxD)	01712	1	299,00	8
SmartRack® starter set Biology		01507	1	1.099,00	11
SmartRack® starter set Chemistry		01508	1	1.099,00	12

## SmartRack® | Accessories

Item	Type	Item no.	PU	Price in EUR	Page
Drip tray	for SmartRack®	01603	1	99,99	19
Drip tray mini	for SmartRack® mini	01604	1	89,99	19
Slot rail mini with handles	for SmartRack® mini	01602	1	54,99	19
Slot rail with handles	for SmartRack®	01601	1	74,99	19
Back-to-back connectors	for SmartRack®	01614	1	79,99	20
Protective cover for SmartRack®	with zipper	01609	1	139,99	21
Protective cover for SmartRack®	without zipper	01613	1	89,99	21
Protective cover for SmartRack® mini	with zipper	01608	1	99,99	21
Protective cover for SmartRack® mini	without zipper	01612	1	64,99	21

## SmartCase

Item	Type	Item no.	PU	Price in EUR	Page
SmartCase		02101	1	949,00	22

Net prices plus statutory VAT. | Valid from 01.01.2026  
 The general terms and conditions of Better Basics Laborbedarf GmbH apply.

# Technical Information

## Technical information:

- All products are washable with water and mild detergent.
- In addition all components can be wiped with alcohol-based disinfectants.
- The plastics and metals used are 100% recyclable.
- Our products are autoclavable up to 134 °C and 3 bar.
- The plastic elements are resistant to heat deformation under normal laboratory conditions.
- To gently clean the anodised aluminium rail, we recommend using a pH-neutral hand-washing detergent. Ensure that the hand wash detergent is odourless and contains no perfumes, greases or additives. Clean the slot rail by hand and rinse thoroughly with clean water.

## Note on joint tolerances for ST14:

- Maximum outside diameter at the narrowest point of the cone joint: 17.2 mm (larger diameters are too wide for the holder and cannot be held)
- Minimum outer diameter at the widest point of the cone joint: 18.0 mm (smaller diameters are too narrow for the holder and would fall out)

## Note on joint tolerances for ST29:

- Maximum outside diameter at the narrowest point of the cone joint: 31.4 mm (larger diameters are too wide for the holder and cannot be held)
- Minimum outer diameter at the widest point of the cone joint: 32.8 mm (smaller diameters are too narrow for the holder and would fall out)

**IMPORTANT TO NOTE:** The products of the SmartIntegrate line, the SmartRack® and SmartRack® mini as well as the SmartCase and SmartStand are not suitable for the permanent storage of hazardous substances.

We reserve the right to make technical changes, material or model changes, errors and misprints excepted. If you have any further questions, please do not hesitate to contact us at [contact@better-basics.de](mailto:contact@better-basics.de).

## For more detailed information, please visit our website:

Functionality of SLS 3D printing:	<a href="http://www.sls-3d-printing.better-basics-laborbedarf.de">www.sls-3d-printing.better-basics-laborbedarf.de</a>
Resistance to UV light:	<a href="http://www.uv-resistance.better-basics-laborbedarf.de">www.uv-resistance.better-basics-laborbedarf.de</a>
Resistance to aggressive media:	<a href="http://www.material-resistance.better-basics-laborbedarf.de">www.material-resistance.better-basics-laborbedarf.de</a>
Sustainability and design-efficient manufacturing:	<a href="http://www.sustainability.better-basics-laborbedarf.de">www.sustainability.better-basics-laborbedarf.de</a>



Photo: Building of the Better Basics Laborbedarf GmbH in Dresden

## YOU ARE INTERESTED IN OUR PRODUCTS OR YOU HAVE FURTHER QUESTIONS?

Get in touch with us:

[contact@better-basics.de](mailto:contact@better-basics.de)

You can also find more information and news about our product innovations at

[www.Better-Basics-Laborbedarf.de](http://www.Better-Basics-Laborbedarf.de)

### **Better Basics Laborbedarf GmbH**

Loebtauer Str. 69

01159 Dresden

Germany

T: +49 (0) 178 9617 577 (Direct customers)

T: +49 (0) 176 6233 8026 (Sales partners)

E: [contact@better-basics.de](mailto:contact@better-basics.de)

W: [www.Better-Basics-Laborbedarf.de](http://www.Better-Basics-Laborbedarf.de)

### **Our request to you**

Our products such as our SmartRack® and our SmartIntegrate products are only as good as their users. With our roots in the Leibniz Institute for Polymer Research Dresden, our company Better Basics Laborbedarf stands for the tradition of “Made in Germany” quality. We feel in multiple aspects committed to this quality label for German companies, which stands for high innovative strength. Even beyond our product catalogue, the same applies for us at Better Basics: Just contact us if you have ideas for extensions or improvements or if you notice something in your daily laboratory routine that we should optimise right away.

We want to create the best possible working environment for you in the laboratory and we are grateful for any comments and inspiration. We can jointly shape the future of laboratory work by cooperating with scientists and researchers. Just like you, we’re always looking for a way to make our world a little better every day.

### **Impress:**

Better Basics Laborbedarf GmbH  
Loebtauer Str. 69, G-01159 Dresden

E-Mail: [info@better-basics.de](mailto:info@better-basics.de)

CEO: Mario Schneider

Commercial register: HRB 39524; Register court: Amtsgericht Dresden

VAT registration number: DE327508812

